



User Manual

Dil-Membrane EVs Labeling & Purification Kit (orange-red)

Cat. # EXOPDil10-1
EXOPDil20-1

辽宁润基生物科技有限公司

Liaoning Rengen Biosciences Co.,Ltd

Contents

Storage and Application.....	2
Product Description.....	3
List of Components.....	3
General Information.....	4
Protocols	
Staining Protocol.....	4
Purification Protocol.....	5
Example Data and Applications.....	6
Related Products.....	7
Troubleshooting	7
Technical Support.....	8

Storage and Application

【Storage】

The Dil-Membrane EVs Labeling & Purification Kit (red) is shipped on ice, the components should be stored at recommended temperatures and protected from light. Properly stored kits are stable for 6 months. please read the instructions before use.

【Application】

For research use only. Not for use in diagnostic procedures.

Product Description

The carbocyanine dye Dil poses highly lipophilic nature and is widely used to stain cell membranes. The dye uniformly labels cells via lateral diffusion in the plasma membrane. The fluorescence of free Dil is very weak, but greatly enhanced when incorporated into membranes. Dil emits orange-red fluorescence when excited, its maximum excitation and emission wavelength is **549nm** and **565nm**, respectively.

Extracellular vesicles (EVs) are membrane-derived particles surrounded by a phospholipid bilayer that are released by cells. We use Dil dye to efficiently label EVs membranes. Then the excess unbound dyes are fast removed through our excellent Spin Columns from labeled exosome preparations, which separate molecules on the basis of the differences in size. Contaminant removal is easier and faster than traditional clean-up methods such as ultracentrifugation, spin filters.

With a highly specific membrane labeling and very low background levels, the kit can be used for most applications that require visualization of labeled EVs for tracking studies.

List of Components

Specification: **20 reactions/Kit (Cat.# EXOPDil20-1);**
10 reactions/Kit (Cat.# EXOPDil10-1)

Item (Cat.#EXOPDil20-1)	Volume	Storage Temp.
Dil Labeling Dye*	100µL	-20°C
Reaction Buffer	1.0mL	2-8°C
Spin Columns# within Collections Tubes	20 sets	2-8°C
1.5mL Light-proof Microcentrifuge Tubes	20 tubes	RT
Item (Cat.#EXOPDil10-1)	Volume	Storage Temp.
Dil Labeling Dye*	50µL	-20°C
Reaction Buffer	500µL	2-8°C
Spin Columns# within Collections Tubes	10 sets	2-8°C
1.5mL Light-proof Microcentrifuge Tubes	10 tubes	RT

Note:

1. * Protect labeling dye from light.
2. # Keep the Spin Columns upright stand on end.
3. The sterilized PBS buffer is not provided in the Kit. Please prepare at least 600uL PBS buffer for each reaction.

General Information

1. This Kit can label exosomes of any resources, including cell culture supernatants and body fluids (such as serum, plasma, urine, CSF or saliva).
2. When extracting exosomes for Labeling, the minimum dosage of serum and plasma is 500μL, the urine is 10mL, and the cell culture supernatant is 5mL.
3. It is not recommended to use exosomes extracted by PEG precipitation method, which contain too many impurities. The exosomes extracted by ultracentrifugation, affinity method or our company's **Exosome Extraction and Purification Kit** (for serum/plasma) or **Exosome Concentration Kit** (for Cell media/urine) is preferable.
4. Heat the dye at 37 °C until dissolved completely when the dye has crystallized.
5. Fluorescent dyes have quenching problems, please protect dyes from light during operating.
6. For your safety and health, please wear a lab gown and wear a disposable glove when operating.

Protocols

Staining Protocol

1. Vortex the DiO Labeling Dye and then instantaneous centrifugation, add 5μL DiO Labeling Dye to 50μL Reaction Buffer, and mix well until the dye is dissolved completely.
2. Add 50μL exosomes samples into dye mixture from step 1, and mix well.
3. Incubate the mixture for 30 minutes at 37 °C, pipet twice during the incubation period.

Note: *Protect the tubes from light.*

Purification Protocol (remove excess unbound dye)

4. Prepare the Spin Column prior to application of your sample.
 - a) Open the lip of Spin Column within Collection Tubes, using a micropipette, aspirate preservative buffer from the top of the column and discard it. In order to prevent drying of the column bed, proceed to the next step promptly.
 - b) Remove the outlet plug, equilibrate the column by adding 200 µL sterilized PBS buffer (not provided) and spin down at 50 x g for 90 seconds. If any PBS remains above the top filters, repeat spin at the same speed with 10 seconds increments, discard the eluate.
 - c) Repeat the procedure b) again.


Note: *Do not spin at too high speed or for too long as this may desiccate or compress the resin and decrease the function of spin column.*

5. Carefully apply 100 µL exosome labeling preparation (from step 3 above) to the top of the column.

Note: *The maximum capacity of the spin column is 100µL. Do not load samples more than 100µL.*

6. Centrifuge at 50 x g for 90 seconds, discard the eluate.
7. Place the column into a 1.5 ml Light-proof Microcentrifuge Tubes (provided). Apply 200 µL PBS buffer (not provided) to the top of the column.
8. Centrifuge at 50 x g for 90 seconds. The 200 µL eluate contains the labeled exosomes.
9. Labeled exosomes can be added to the cells or used for downstream applications.

Example Data and Applications

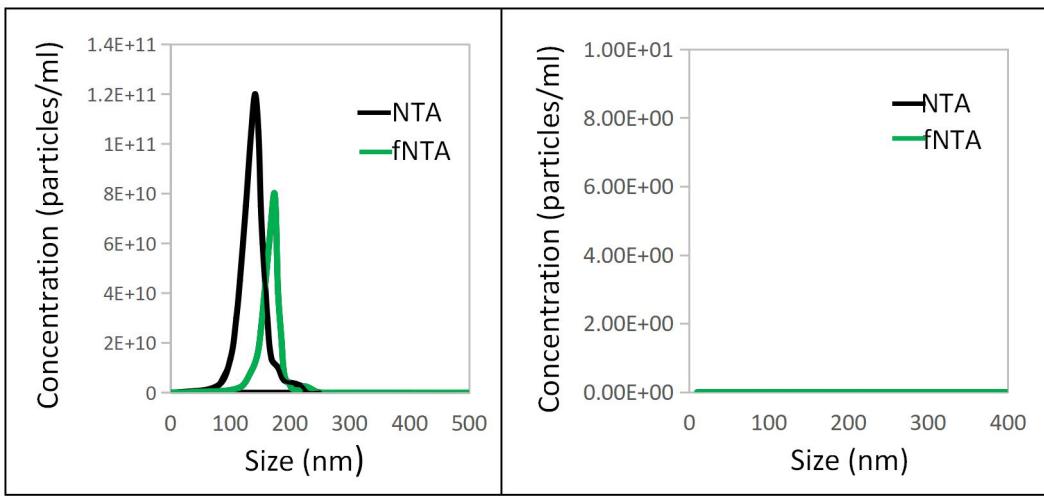

A: Dil labeled EVs

B: Dye only (control)

Figure 1. Dil labeled EVs and free dye control were analysed by NTA and fNTA.

The exosomes from LnCAP cell culture supernatant were concentrated by Exosome Concentration Kit (Cat# EXOCCon05-10), then were labeled and purified with Dil Membrane EVs Labeling & Purification Kit. A control involving dye but no exosomes was performed in parallel to confirm dye retention by the column. The results of NTA and fNTA indicated that a fairly considerable portion of exosomes were successfully labeled, the free dyes did not aggregate and the unbound dye can be removed successfully through the column.

Related Products

Exosome labeling & Purification	
DiO-Membrane EVs Labeling & Purification Kit (green)	EXODiO10-1/EXODiO20-1
Dil-Membrane EVs Labeling & Purification Kit (orange-red)	EXODil10-1/EXOIl20-1
DiR-Membrane EVs Labeling & Purification Kit (near-infrared red)	EXODiR10-1/EXOIr20-1
PKH67-Membrane EVs Labeling & Purification Kit (green)	EXOPKH67-10/EXOPKH67-20
Exosome Isolation & Purification	
Exosome Extraction & Purification Kits (for blood serum/plasma)	EXORG10SP-1/ EXORG30SP-1/
Exosome Concentration Kits (for cell culture media/urine)	EXOCCon5-10/ EXOUCon5-10
Total Exosome Capture & Isolation Kits (for cell culture media/urine)	EXOMBoCU-10/EXOMBoCU-20
Exosome Nucleic Acid Extraction	
Exosome Extraction & DNA Isolation Kits (for blood serum/plasma)	EXODNA30A-1/ EXODNA50A-1
Exosome Extraction & DNA Isolation Kits (for cell culture media/urine)	EXODNA10B-1/EXODNA24B-1
Exosome Extraction & RNA Isolation Kits (for blood serum/plasma)	EXORNA30A-1/EXORNA50A-1
Exosome Extraction & RNA Isolation Kits (for cell culture media/urine)	EXORNA10B-1/EXORNA24B-1
Exosome Spin Columns (MW4000)	RGESC12-1/RGESC24-1

Troubleshooting

Q1: The fluorescence signal is low than expected?

A1: There may be too low amount exosome for labeling, it is recommended to take a larger sample to extract exosome. In addition, fluorescent dye will be quenched, please keep dye from light when operating. Ensure that the columns do not dry out during the procedure. Spinning the column for too long or at too high speed may cause the column to work inefficiently and loss of the labeled exosomes.

Q2 : How to store the labeled exosomes when not carry out the downstream experiments immediately?

A2: The labeled exosomes can be stored at 2-8°C for 1-2 days aware of light, If kept for a long time, it is recommended to store at -80 °C avoid repeated freezing and thawing.

Q3: Can I increase the elution volume?

A3: This is not recommended as it will result in co-elution of excess unbound dye with exosomes.

Q4: How I can know the counts of labeled exosomes?

A4 : The labeled exosomes usually can be detected and counted by fluorescent Nanoparticle Tracking Analysis (fNTA) or flow cytometry. Dil labeling did not change the size of exosomes actually, but the calculation of particle size is related to many factors, which leads to the measurement value of size of the same labeled exosome sample under the fluorescence mode (fNTA) is larger than that under the conventional mode (NTA).

Technical Support

For more information about our products and to download manuals, please visit our web site: <http://www.rengenbio.com>

For additional information or technical assistance, please call or email us.

Liaoning Rengen Biosciences Co., Ltd.

Add.: Building 20th, LIANDO U Valley, Number 77 Road 13th, Shenyang Economic and Technological Development Zone, Liaoning Province, 110027

Phone: 024-3108,6590

Fax : 024-3108,6589

E-mail: General Information: info@rengenbio.com

Technical Support: support@rengenbio.com

Ordering Information: order@rengenbio.com



Wechat Public Platform



Exosome Research Exchange Group